

Food Allergens

Veratox®

For Total Milk Allergen



Intended Use

Veratox for Total Milk Allergen is used for the quantitative analysis of milk residue in food products such as juices, wine, sauces or sorbets.

The Test

The test is a sandwich ELISA that provides exact concentrations in parts per million (ppm). Any present milk residue is extracted from samples with a buffered salt solution (PBS) by shaking in a heated water bath. Extracted milk residue is sampled and added to antibody-coated wells (capture antibody). Any unbound residue is washed away and a second, enzyme-labeled antibody (detector antibody) is added. The detector antibody binds to the already bound milk residue. After a second wash, the substrate is added. Color develops as a result of the presence of bound detector antibody. Red Stop reagent is added and the color of the resulting solution is observed. A microwell reader is used to yield optical densities. Control optical densities form a standard curve, and sample optical densities are plotted against the curve to calculate the exact concentration of milk residue.


The Procedure

Samples must be extracted prior to testing.

1. Add 150 μ L controls and extracted samples to transfer wells.
2. Transfer 100 μ L to the antibody wells. Incubate for 10 minutes.
3. Dump liquid from antibody wells.
4. Wash wells thoroughly with wash buffer. Tap out water on paper towel.
5. Transfer 100 μ L conjugate from reagent boat to antibody wells using 12-channel pipettor. Incubate for 10 minutes.
7. Repeat steps 3-4 by dumping out the liquid, thoroughly washing the wells, and tapping dry.
8. Transfer 100 μ L substrate from reagent boat to antibody wells using 12-channel pipettor. Incubate for 10 minutes.
9. Transfer 100 μ L Red Stop from reagent boat to antibody wells.
10. Read results in a microwell reader.

See package insert for complete instructions

Results

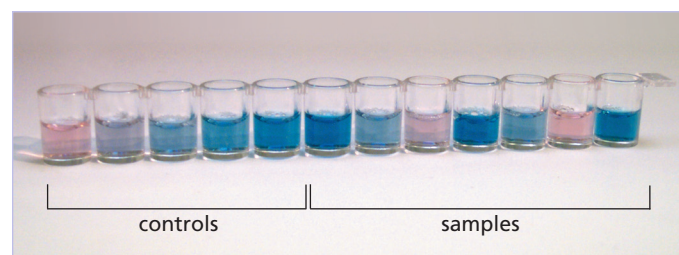

Product Specifications

Range of quantitation:	2.5 ppm - 25 ppm
Controls provided:	0, 2.5, 5, 10 and 25 ppm non-fat dry milk
Testing time:	30 minutes
Tests per kit:	Up to 38

Materials Recommended But Not Provided

Available from Neogen*

1. Allergen Extraction Kit
2. Adjustable 50-200 μ L pipettor
3. Graduated cylinder capable of measuring 125 mL
4. Scale capable of weighing 5 ± 0.1 g
5. Microwell reader with a 650 nm filter
6. 50-200 μ L pipettor
7. 12-channel pipettor
8. Tips for adjustable and 12-channel pipettors
9. Timer
10. Microwell holder
11. Wash bottle
12. 3 reagent boats for use with 12-channel pipettor
13. 1.0 L bottle to prepare washing solution
14. 1.0 L heat safe bottle to prepare extract solution

Not available from Neogen

1. Shaker water bath adjusted to 60°C (140°F) with clamps to hold extraction bottles
2. Paper towels or equivalent absorbent material
3. Waterproof marker
4. Distilled or deionized water

Ordering Information

Prod.#	Product description
8470	Veratox for Total Milk Allergen
8429	Allergen Extraction Kit - 20 samples